

Fixvo II Microlitre Pipette

**LA599, LA600, LA601, LA602,
LA603, LA604, LA605, LA606,
LA607, LA608, LA609, LA610**

These are fixed general purpose micropipettes for sampling and dispensing accurate amounts of liquid in micro volume. Pipettes operate on the air displacement principle and use disposable tips. All models mentioned below are equipped with a built-in tip ejector.

A range of economic, accurate, fixed volume and one-hand dispensing of fixed volume in routine laboratory procedures micro-pipettes.

Application : Ideal for molecular biological applications, Medical laboratories, Medical research and industrial laboratories etc.

Cleaning of Pipette: To clean the pipette use ethanol and a soft cloth or lint-free tissue. It is recommended to clean the tip cone regularly.

Calibration : Each pipette has been checked and calibrated at factory with procedure conforming to DIN 12650 standards. It is recommended to check the calibration atleast once a year , for regularly used pipette.

Storage : When not in use it is recommended that the pipette be stored in a vertical position. Keeping pipette on its side can cause liquids to leak in to the body of the pipette and cause corrosion.

Product Name	Product Code	Description	Capacity (µl)
Fixvo II Micropipette-5	LA599	Fixvo II Micropipette	5.00
Fixvo II Micropipette-10	LA600	Fixvo II Micropipette	10.00
Fixvo II Micropipette-20	LA601	Fixvo II Micropipette	20.00
Fixvo II Micropipette-25	LA602	Fixvo II Micropipette	25.00
Fixvo II Micropipette-50	LA603	Fixvo II Micropipette	50.00
Fixvo II Micropipette-100	LA604	Fixvo II Micropipette	100.00
Fixvo II Micropipette-200	LA605	Fixvo II Micropipette	200.00
Fixvo II Micropipette-250	LA606	Fixvo II Micropipette	250.00
Fixvo II Micropipette-500	LA607	Fixvo II Micropipette	500.00
Fixvo II Micropipette-1000	LA608	Fixvo II Micropipette	1000.00
Fixvo II Micropipette-2000	LA609	Fixvo II Micropipette	2000.00
Fixvo II Micropipette -5000	LA610	Fixvo II Micropipette	5000.00

Disclaimer :



User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.