

MBT088

ChromTaq DNA Polymerase

Components

Reagents provided	MBT088		
	100 Units	200 Units	500 Units
ChromTaq DNA Polymerase (1 U/μl)	100 μl	200 μl	500 μl
10X HiBuffer A (without MgCl ₂)	400 μl	800 μl	2 ml
10X HiBuffer S (with 17.5 mM MgCl ₂)	400 μl	800 μl	2 ml
50mM MgCl ₂	200 μl	400 μl	1 ml

Description:

Taq DNA Polymerase is a thermostable DNA polymerase of thermophilic bacterium *Thermus aquaticus*. It is suitable for applications requiring high temperature synthesis of DNA. The enzyme catalyzes 5'-3' synthesis of DNA, has no detectable 3'-5' exonuclease (proof-reading) activity, but maintains the 5'-3' exonuclease activity. In addition, Taq DNA Polymerase exhibits deoxynucleotidyl transferase activity, resulting in addition of extra adenines at 3'-end of PCR products.

Features:

- Color tracer dyes for ease of visualization of the addition of polymerase to the reaction.
- Use of 1 U in every 25 μl final reaction allows direct loading and serves as tracking dyes during gel electrophoresis. The blue and pink color dyes migrate approximately at 4 kb and 0.3 kb respectively on 1% TAE agarose gel.
- Thermostable enzyme of approximately 94 kDa from *Thermus aquaticus*.
- Ultra pure recombinant protein.
- Replicates DNA at 74°C and exhibits a half-life of 40 minutes at 95°C.
- Generates mostly 3' dA overhang PCR products which are suitable for TA cloning.
- DNA sequencing and labeling.
- Incorporates modified nucleotides (E.g.: biotin, deoxygenin, fluorescently-labelled nucleotides).
- Supplied with two buffers- **10X Hibuffer A** with KCl and **10X Hibuffer S** with (NH₄)₂SO₄. [(NH₄)₂SO₄ allows for PCR at wide range of Mg²⁺ concentrations and decrease unspecific priming].

Concentration: 1 U/μl

Source: *E.coli* cells with a pol gene from *Thermus aquaticus*

Molecular weight: 94 kDa monomer

Unit Definition:

1U is defined as amount of enzyme that is required to catalyze the incorporation of 10 nmoles of dNTP into acid-insoluble material in 30 minutes at 74°C.

Reaction Buffer:**10X HiBuffer A (Without MgCl₂)**

500mM KCl, 100mM Tris-HCl (pH 9.1 at 20°C) and 0.1% Triton X-100. The buffer is optimized for use with 0.1-0.2mM of each dNTP.

10X HiBuffer S

160mM (NH₄)₂SO₄, 500mM Tris-HCl (pH 9.2 at 22°C), 17.5mM MgCl₂ and 0.1% Triton X-100. The buffer is optimized for use with 0.35mM of each dNTP.

Guidelines for PCR optimization using HiMedia's ChromTaq DNA Polymerase:

- **DNA Template**
 1. Use high quality, purified DNA templates.
 2. Approximately, 10⁴ copies are required to detect the amplification in 25-30 PCR cycles.
 3. Use higher DNA concentration when few PCR cycles are desired.
- **Primers**
 1. Generally 20-30 bp in size.
 2. GC content between 40-60% ideally.
 3. Melting temperatures should be between 42-65°C.
 4. Final concentration to be used 0.1-0.5µM of each primer.
- **Magnesium Concentration**
 1. Ideal for ChromTaq DNA Polymerase is 1.5-2.0mM.
 2. Optimum concentration depends on template, buffer and dNTPs.
 3. Higher than optimal concentration yields undesired products and if concentration is too low the concentration, no amplification products are detected.
- **dNTPs**
 1. Typical concentration to be used is 200µM.
 2. Higher than optimal concentration of dNTPs yields higher yield but low fidelity.
- **ChromTaq DNA Polymerase**

Typical concentration to be used is 0.5 to 2 units per 50µl of reaction.
- **PCR reaction**
 1. Thaw all reaction components on ice.
 2. To PCR reaction, add Taq Polymerase at the end.
 3. Once the reaction is set, immediately transfer the tubes to pre-heated thermal cycler.
 4. Start the reaction with desired cycling conditions with annealing temperature set to 5°C difference of melting temperature between forward and reverse primers.

Quality Control:

All preparations are assayed for contaminating endonuclease, exonuclease and non-specific DNase activities. Functionally tested in DNA amplification.

Buffers recommended for different sizes of template DNA

Buffers	Size of template DNA to be amplified		
	100bp-5kb	5kb-8kb	8kb-20kb
HiBuffer S (1X)	-	+	+
HiBuffer A (1X)	+	-	-
MgCl ₂	+	+	+

Key: + Indicates recommended buffer

Inhibition and Inactivation:

- Inhibitors: ionic detergents (deoxycholate, sarkosyl and SDS) at concentrations higher than 0.06, 0.02, and 0.01% respectively.
- Inactivated by phenol/chloroform extraction.

NOTE:

The error rate for ChromTaq DNA Polymerase, which lacks proof-reading activity is approximately 1 to 2 X 10⁻⁵ errors (or mutation frequency) per nucleotide per duplication. Accordingly, the accuracy of PCR is 4.5X10⁴. Accuracy is an inverse of the error rate and shows an average number of correct nucleotides incorporated before an error occurs.

Storage conditions: The ChromTaq DNA Polymerase should be stored at -20°C. When stored under the recommended conditions, the product is stable for 2 years.

Representative data:

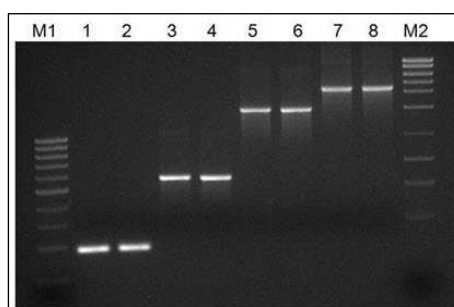


Figure representing amplification of different amplicon sizes using ChromTaq DNA Polymerase with HiBuffer A and HiBuffer S.

Technical Assistance

At HiMedia's, we pride ourselves on the quality and availability of our technical support. For any kind of technical assistance, mail at mb@himedialabs.com

PIMBT088_0/0714 MBT088-01

Disclaimer :

User must ensure of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ Publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.