

Bradford Kit for Protein Estimation

<u>Product Name</u>	<u>Product Code</u>	<u>Kit Packing</u>
Bradford Kit for Protein Estimation	ML178-1PK	1pk

Introduction: HiMedia's Bradford Kit for Protein Estimation is formulated for rapid and accurate quantitative estimation of protein samples by Bradford assay method which was developed by Marion M. Bradford in 1976. This kit contain ready-to-use Bradford Reagent and Protein Standard.

Description: Bradford Reagent contains a dye, Coomassie brilliant Blue G-250, which has an absorbance maximum of 465 nm in unbound state. The Bradford protein assay is based upon the formation of complexes between Coomassie Brilliant Blue G-250 and the protein samples in solution. When the protein sample binds to the dye the color of the solution turns blue from brown and there is a shift in the absorption maximum of the dye from 465 nm to 595 nm. This dye binding procedure is completed within 5 minutes and the blue colored complex formed is stable for one hour. Thus concentration of unknown protein sample can be derived by plotting its absorbance value on the standard curve. The standard curve is obtained from the absorbance readings of the series of standard protein dilutions assayed alongside the unknown sample.

Application: HiMedia's Bradford Reagent provide accurate determination of protein concentration with most of the protein samples. Following are the features of Bradford Reagent.

1. Samples can be estimated visually or values can be determined with a standard Spectrophotometer or plate reader at 595nm
2. HiMedia's Bradford reagent is ready-to-use and no need to prepare working standards.
3. Bradford assay is much easier and faster than traditional Lowry method.
4. Detects protein concentration in the range 20 to 1500 µg/mL

Kit Contents

Product Code	Reagents	Quantity	Storage
ML106	Bradford Reagent	500 ml	2 – 8°C
MBT129	Protein Standard (1 mg/ml)	1X5 ml	2 – 8°C

Materials required but not provided: Spectrophotometer, Test tubes, 1 ml disposable cuvettes, 96-well plate.

Storage conditions: Bradford Reagent and Protein Standard should be stored at 2 - 8 °C

Estimation of unknown protein using standard curve (with BSA)

1. Take seven test tubes and label them as Blank and 1 to 6.
2. Make dilutions of standard protein (BSA) with concentrations of 20, 16, 12, 8, 4 $\mu\text{g}/200 \mu\text{l}$ by transferring respective amount of BSA solution (stock: 1mg/ml) and adjusting it to a total volume of 200 μl by adding distilled water as mentioned in the following table.
3. Add 1ml of Bradford's Reagent to each test tube and mix the contents of each tube thoroughly by vortexing the tubes and incubate at room temperature (below 30°C) for 10 minutes.
4. Transfer the content of the tubes to cuvettes and measure the absorbance at 595 nm wavelength.

Tube No.	Blank	1	2	3	4	5	
Conc. of BSA (μg)	0.0	4	8	12	16	20	Unknown sample
Amt of Stock (μl)	0.0	4	8	12	16	20	200 μl
Amt of diluent (μl)	200	196	192	88	84	80	
Amt of Bradford Reagent (ml)	1	1	1	1	1	1	
Vortex the tubes and incubate at RT for 10 minutes							
Absorbance at 595 nm							

5. Plot a Standard Curve of absorbance at 595 nm on "Y" axis versus concentration of protein $\mu\text{g}/200 \mu\text{l}$ on "X" axis.
6. Record the value "x" of unknown sample from graph corresponding to the absorbance reading.

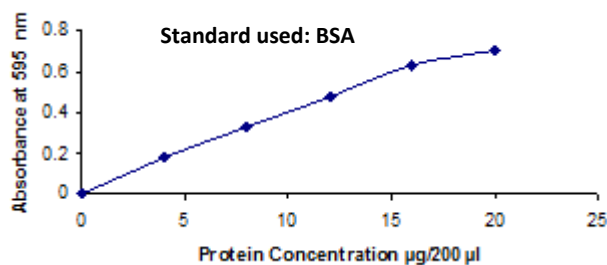


Fig: Standard curve for protein estimation by Bradford assay

Protein concentration can be calculated using following formula:

$$\text{Protein Concentration in Unknown Sample} = \frac{\text{Concentration of Unknown in } \mu\text{g}}{\text{Volume of sample in } \mu\text{l}} \times 1000 \mu\text{g/ml}$$

Please refer disclaimer Overleaf.

96-well plate assay:

This assay should be performed in a 96-well plate. It is very fast and easy to quantitate multiple protein samples at a time.

NOTE: For Standard Graph, 5 μ l of 0.2-1.4mg/ml of protein sample can be used.

Protocol:

1. Prepare protein standards as mentioned above
2. Add 5 μ l of standard protein samples to each well of 96-well plate.
3. Add 5 μ l of unknown samples to the wells.
4. To each well, add 250 μ l of Bradford reagent.
5. Mix well by pipetting or keep it in a shaker for 15-30 seconds.
6. Incubate the samples for 5-45 minutes at room temperature and measure the absorbance at 595nm.
7. The absorbance readings must be taken before 60 minutes or within 10 minutes after the completion of incubation.
8. Plot a Standard Curve of absorbance at 595 nm on “Y” axis versus concentration of protein on “X” axis.
9. Record the value “x” of unknown sample from graph corresponding to the absorbance reading and determine the protein concentration.

Technical Assistance:

At HiMedia we pride ourselves on the quality and availability of our technical support. For any kind of technical assistance, mail at mb@himedialabs.com.



Consult instructions for use



Do not use if package is damaged



HiMedia Laboratories Pvt. Limited,
Reg. Off: 23 Vadhani Industrial Estate,
LBS Marg, Mumbai - 400086,
India Works: B/4-6, M.I.D.C., Dindori,
Nashik, India (or respective plant address)
Customer Care No: 022-6116 9797
www.himedialabs.com

PIML178_0/07 17

ML178-00

Disclaimer :

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.

HiMedia Laboratories Pvt. Ltd. A-516, Swastik Disha Business Park, Via Vadhani Ind. Est., LBS Marg, Mumbai-400086, India.
Customer care No.: 022-6147 1919 Email: techhelp@himedialabs.com Website: www.himedialabs.com