

## HiDispo™ Bag

**PW038, PW040, PW048,  
PW055, PW056 & PW057.**

These bags are autoclavable bag used for waste disposal and recommended for disposal of pathological/ clinical or contaminated material. It can also be used for sterilization of glass ware or plastic wares.

**Application:** Used in hospitals, clinical laboratories, microbiological laboratories, Biotechnology laboratories, molecular biology laboratories etc. for disposal of clinical and pathological or contaminated material.

### Direction for use:

1. Place all the clinical / contaminated material in the HiDispo bag.
2. Seal the bag.
3. Autoclave at 121°C for 15 minutes.
4. Discard the autoclaved HiDispo bag with decontaminated material as per local law.

Product Name	Product Code	Description	Size (Inch) H X B	Packing
HiDispo™ Bag-14	PW038	Clear, transparent and autoclavable disposable bag.	20" X 14"	500 & 1000 Nos.
HiDispo™ Bag-12	PW040	Clear, transparent and autoclavable disposable bag.	12" X 10"	500 & 1000 Nos.
HiDispo™ Bag-10	PW048	Clear, transparent and autoclavable disposable bag.	10" X 6"	500 & 1000 Nos.
HiDispo™ Bag-40	PW055	Clear, transparent and autoclavable disposable bag.	40" X 20"	250 & 500 Nos.
HiDispo™ Bag-34	PW056	Clear, transparent and autoclavable disposable bag.	34" X 20"	250 & 500 Nos.
HiDispo™ Bag-7	PW057	Clear, transparent and autoclavable disposable bag.	11" X 7"	500 & 1000 No

### Products features:

- Superior strength polypropylene bag, are free from chlorine.
- Can be steam sterilization at 121 °C- 15 minutes.
- Decontamination and Disposal of contaminated material made easy.
- Transparent – easy to see through.
- Available in multiple packing of 50 bags.